

Micro Distillations: SimpleDist® Micro Distillation Block Environmental Express® Disposable Distillation Micro Tubes


Eliminate costly glassware and increase the speed of distillations in your lab!

- Distill up to 26 samples simultaneously
- Collection tubes snap easily at score line
- Molded-in graduations at 6mL and 10mL make measuring easy

Each box of Environmental Express disposable distillation micro tubes contains:

- A. Collection tubes (shown w/cap)
- B. Corresponding number of sample collection tubes
- C. Caps
- D. Hydrophobic membranes
- E. Complete tube assembly.

A variety of pack sizes are available to meet the needs of your laboratory.

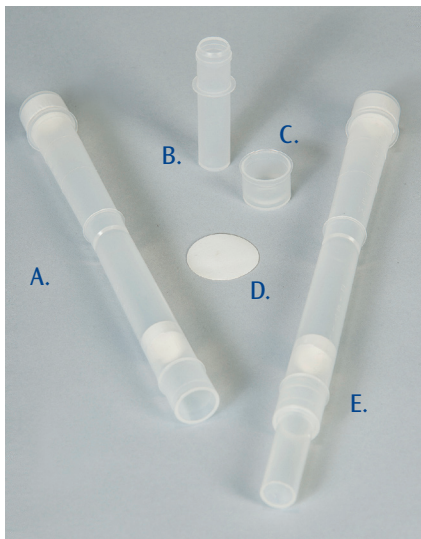

Environmental Express Disposable Distillation Micro Tubes Instructions for Use:



1. Pipet 6mL of sample into each sample tube.
2. Securely attach the sample tube to the collection tube.
3. Place the assembled tubes into the SimpleDist Micro (or other micro distillation block) and heat the samples according to method.



4. Wearing heat protective gloves, carefully separate the sample tube from the collection tube and cool in rack.



5. Snap the collection tube in two at the score line.



6. Bring contents of collection tube up to volume with DI Water.

**Economical Tubes
Fit Your Existing
Microdistillation Block!**

Please download your SimpleDist Micro Product Manual from the Technical Support/Product Manuals section of our website prior to operating your block

ENVIRONMENTAL EXPRESS

Call 800.343.5319 or 843.881.6560 • www.environmentalexpress.com

